
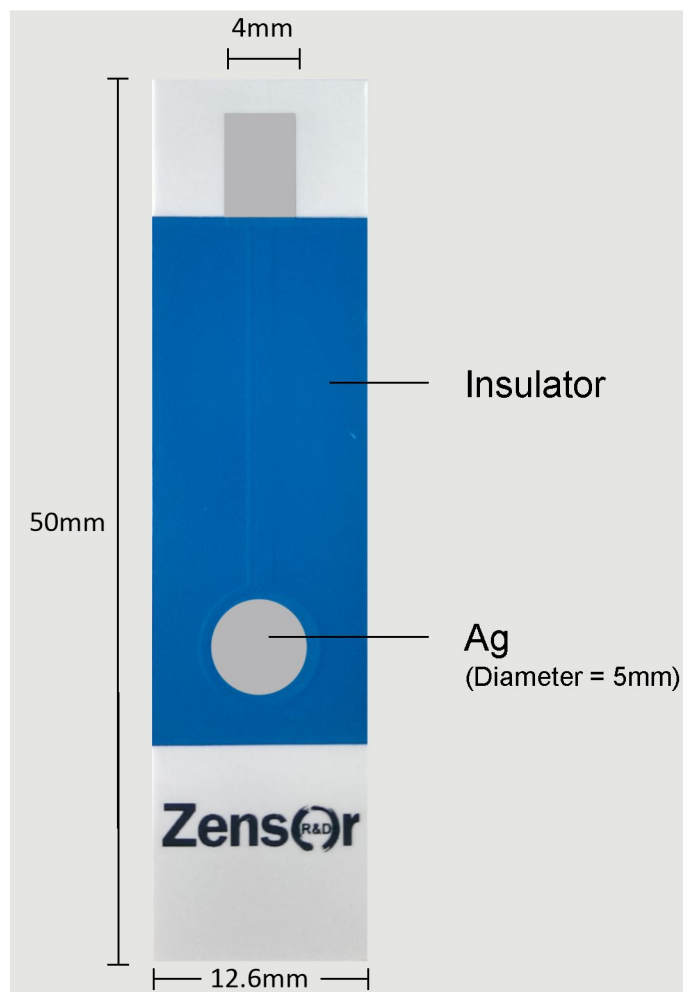


### PRODUCT DESCRIPTION

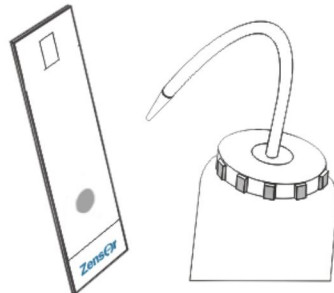

### Before Use (Recommend)

- A** Flush it with DI water for several times.
- 
- B** Immerse it in DI water for 5~10 min.
- C** Put it in your buffer solution or 0.1M phosphate buffer solution (pH7.0) and then operation of cyclic voltammetry (CV) with several segments.

### DESCRIPTION FOR USE

1. ASPE is supplied ready for use under acid ( $< 2 \text{ M H}_2\text{SO}_4$ ) and basic media ( $< 1 \text{ M NaOH}$ ).
2. Avoid bending hardly as operating.
3. Should Keeping apart from organic media.

### Storage

Store product in a dry location apart from sunlight at optimal temperature of 20 to 40 °C.