



PRODUCT DESCRIPTION

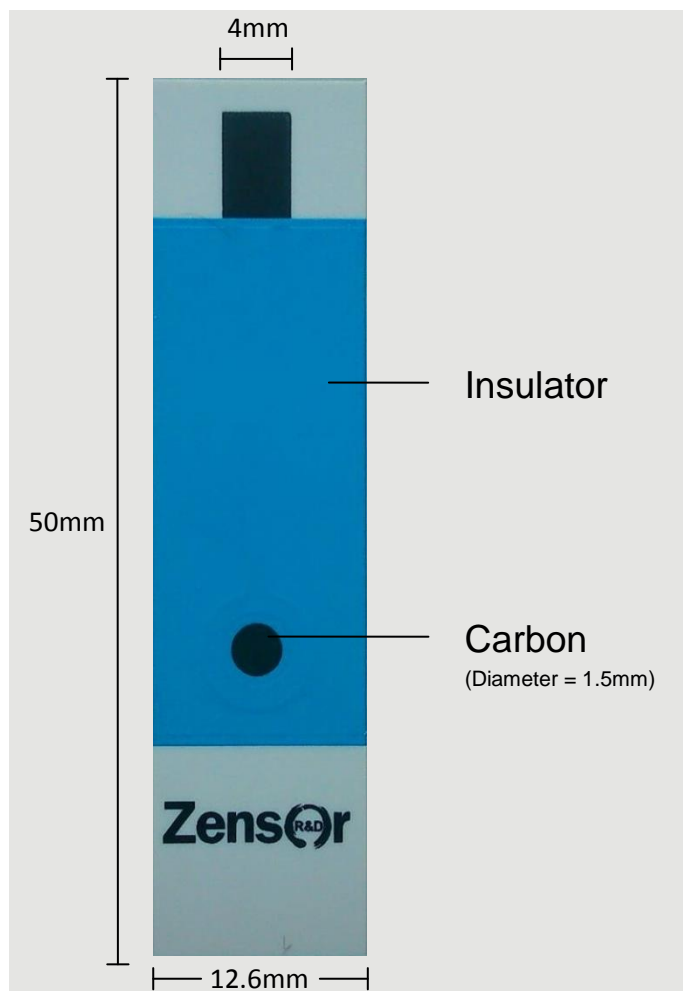

Before Use (Recommend)

A You can wash DI water with several times.

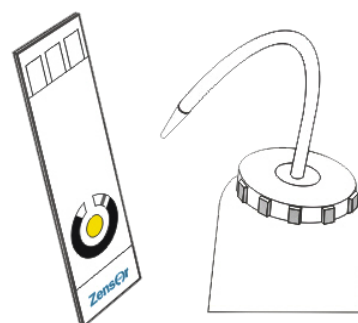

B Or put electrode of SPE into DI water or buffer solution of you prepared at 5~10 min.

C Or put electrode of SPE in 0.1M phosphate buffer solution (pH7.0) and then operation of cyclic voltammetry (CV) with sever segments.

DESCRIPTION FOR USE

1. SE102 is supplied ready for use under acid ($< 2 \text{ M H}_2\text{SO}_4$) and basic media ($< 1 \text{ M NaOH}$).
2. Avoid bending hardly as operating.
3. Should Keeping apart from organic media.

Storage

Store product in a dry location apart from sunlight at optimal temperature of 20 to 40 °C.